



Using deconex[®] products in pipette cleaners

Residue? Inaccurate analytical results?

Avoid this risk by using professional deconex[®] cleaners. Refuse to compromise. deconex[®] products have been specifically developed for cleaning laboratory glassware and ensure residue-free cleaning with consistent and accurate analytical results. deconex[®] cleaning processes offer maximum performance, quality and reliability in the cleaning of laboratory glassware.

Advantages of deconex[®]

Reliable

- Consistent and accurate analytical results
- Excellent rinsability

Professional

- Products developed specifically for laboratory glassware
- Synergetic effect of carefully selected raw materials
- Free from critical ingredients

Economical

- Very cost-effective
- Wide ranging efficacy

deconex[®]


The perfect solutions for research, industry and education.

The following products are recommended for specific applications on the basis of practical tests with Hölzel pipette cleaners in Borer Chemie AG's test laboratories:

Application / dirt type	Product	Dosage	Program
Standard use of pipettes for analysis. Residue from reagents and chemicals. Light to medium soiling	deconex® 15 NF	50 ml	Program 1 40 °C
Proteins, blood Fresh or baked-in dirt Medium to heavy soiling	deconex® 12 PA-x	50 ml	Program 2
Greasy and oily dirt from the pharmaceuticals, cosmetics or petrochemical industry Heavy to very heavy soiling	deconex® FORMULA 1	100 ml	Program 2 90 °C

How can you achieve more process reliability when cleaning pipettes?

What you need to do is qualify and validate your laboratory glassware cleaning process. Let us show you how we can make your work a lot easier.



Help with validation tasks:

- Adopt our validated analysis methods
- Our sample logs make the qualification process easier for you

When qualification and validation are complete, you can be sure that your pipette cleaner, cleaning processes and cleaning agents are effective and optimally suited to each other.

Contact us to find out how much our range can verifiably increase your process reliability.

Borer Chemie AG

Gewerbstrasse 13, 4528 Zuchwil / Switzerland
Tel. +41 32 686 56 00, Fax +41 32 686 56 90
www.borer.ch, office@borer.ch

deconex®